

Physical Developer

Table of Contents

1	INTRODUCTION/SCOPE	2
2	STANDARDS AND CONTROLS	2
3	LIMITATIONS	2
4	EQUIPMENT	2
5	PROCEDURE	2
5.1	Solution Preparation	2
5.1.1	Maleic Acid solution	2
5.1.2	Redox solution.....	2
5.1.3	Detergent solution	3
5.1.4	Silver Nitrate solution.....	3
5.1.5	Physical Developer working solution.....	3
5.1.6	Bleach solution (optional).....	3
5.2	Application	3
5.2.1	Bleach Solution Rinse (Optional)	3
5.3	Storage	4
5.4	Shelf Life.....	4
6	SAFETY	4
7	REVISION HISTORY	4

Physical Developer

1 INTRODUCTION/SCOPE

Physical Developer is used by FBI Laboratory Friction Ridge Discipline personnel to develop latent prints on porous and semi-porous surfaces.

2 STANDARDS AND CONTROLS

See *Processing Overview* ([FRD-300](#)).

3 LIMITATIONS

- A. All metal items, such as staples and paper clips, must be removed from item(s) prior to Physical Developer processing.
- B. Metal tweezers cannot be used during processing.

4 EQUIPMENT

- Citric Acid
- Maleic Acid
- Silver Nitrate
- Bleach
- Synperonic N or Tween 20
- Ferrous Ammonium Sulfate
- n-Dodecylamine Acetate
- Ferric Nitrate

5 PROCEDURE

5.1 Solution Preparation

Personnel will prepare the solutions as follows. Alternative amounts may be prepared, provided the same ratio of chemicals mixed is retained.

5.1.1 Maleic Acid solution

- A. Combine:
 - Maleic Acid - 25 g
 - Distilled water - 1000 mL
- B. Stir until solid dissolves.

5.1.2 Redox solution

In the order listed below, stir until each solid is dissolved before adding the next solid:

- Distilled water - 1000 mL
- Citric Acid - 20 g
- Ferric Nitrate - 30 g
- Ferrous Ammonium Sulfate - 80 g

5.1.3 Detergent solution

- A. Combine:
 - n-Dodecylamine Acetate - 3 g
 - Synperonic N or Tween 20- 4 g
 - Distilled water - 1000 mL
- B. Stir until all chemicals dissolve.

5.1.4 Silver Nitrate solution

- A. Combine:
 - Silver Nitrate - 200 g
 - Distilled water - 1000 mL
- B. Stir until solid dissolves.

5.1.5 Physical Developer working solution

- A. While stirring, combine in the order listed:
 - Redox solution - 1000 mL
 - Detergent solution - 40 mL
 - Silver Nitrate solution - 50 mL
- B. Stir for at least three minutes.

5.1.6 Bleach solution (optional)

Combine:

- Bleach - 500 mL
- Distilled water - 500 mL

5.2 Application

Personnel will complete the following steps in order:

- A. Immerse item(s) in Maleic Acid solution.
- B. Agitate solution, manually or with orbital shaker, for a minimum of 15 minutes.
- C. Immerse item(s) in Physical Developer working solution.
- D. Agitate solution, manually or with orbital shaker, for 10-15 minutes.
- E. Immerse item(s) in first water rinse for at least 1 minute.
- F. Rinse item(s) in second water rinse.
- G. Dry item(s) in air or by applying heat with an iron, heater, or dryer.
- H. Capture appropriate friction ridge detail as applicable (digitally or photographically).

5.2.1 Bleach Solution Rinse (Optional)

- A. The bleach solution may darken latent prints developed with Physical Developer, lighten the background, and remove any previous processing stains that may still be present on the item.
- B. The bleach solution is especially effective on paper bags and paper currency.

- C. The solution is usually applied after the second water rinse but can be done after the item has been dried.
- D. Personnel will complete the following steps in order:
 1. Immerse item(s) in Bleach solution for 10-15 seconds or until desired contrast is achieved.
 2. Rinse item(s) in water.
 3. Dry item(s) as described above.

5.3 Storage

- A. Maleic Acid, Redox, Detergent, and Bleach solutions may all be stored in any type of laboratory acceptable receptacle.
- B. Silver Nitrate solution must be stored in a dark bottle.
- C. Physical Developer working solution is not stored. It is prepared as needed.

5.4 Shelf Life

- A. Maleic Acid, Redox, and Bleach solution have indefinite shelf lives provided the reagent checks are satisfactory.
- B. Detergent and Silver Nitrate solution each have a shelf life of 1 year provided the reagent checks are satisfactory.
- C. Physical Developer working solution is not retained. It is prepared as needed.

6 SAFETY

See [FBI Laboratory Safety Manual](#) for appropriate information.

7 REVISION HISTORY

Revision	Issued	Changes
03	07/15/2021	Replace Latent Print Units with Friction Ridge Discipline. Minor wording changes. Streamline equipment list. Re-organization and re-numbering of sections. Section 4.4 - broken into Section 4.1.1, Section 4.1.2, Section 4.1.3, Section 4.1.4, Section 4.1.5, and Section 4.1.6. Parts of Section 4.2 - broken into Section 4.2.1 and Section 4.2.2. Section 4.1 - added ratio allowance. Section 4.1.6 - moved note. Section 4.2 - removed note. Section 5 - added Preamble
04	07/01/2022	Format Updated. Section 1 – Scope updated.